

# XCAMLITE4K Series Camera Help Manual



## Contents

Contents .....	I
1 The Application of the XCAMLITE4K Series Camera.....	1
2 XCAMLITE4K Series Camera’s Datasheet and Functions(2) .....	1
3 Dimension of XCAMLITE4K Series .....	3
4 XCAMLITE4K Series Camera Packing Information(TBD) .....	3
5 Software and App .....	4
6 XCAMLITE4K Series Camera Application Configurations .....	5
6.1 Camera Working Standalone with Built-in XCamView Software .....	5
6.2 Connecting camera to the PC with USB port.....	5
7 Brief Introduction of XCAMLITE4K Series Camera’s UI and Its Functions .....	6
7.1 XCamView UI .....	6
7.2 The Camera Control Panel on the Left Side of the Video Window .....	7
7.3 The Measurement Toolbar on Top of the Video Window .....	7
7.4 Icons and Functions of the Synthesis Camera Control Toolbar at the Bottom of the Video Window .....	9
8 Sample Photos Captured with XCAMLITE4K Series Camera .....	15
9 Contacting Customer Service .....	15

## 1 The Application of the XCAMLITE4K Series Camera



Figure 1-1 The XCAMLITE4K Series Camera

The XCAMLITE4K series camera is intended to be used for the acquisition of digital images from the stereo microscope and biological microscope. The basic characteristic is listed as below:

- Sony Exmor back illuminated CMOS sensor
- 4K HDMI/USB multiple video outputs
- 4K/1080P auto switching according to the display resolution
- SD card for the captured image and video storage
- Embedded XCamView for the control of the camera
- With strong ISP and other related processing functions
- ToupView/ToupLite software for PC
- ToupLite software for MAC

## 2 XCAMLITE4K Series Camera's Datasheet and Functions(2)

Order Code	Sensor & Size(mm)	Pixel(μm)	G Sensitivity	FPS/Resolution	Binning	Exposure(ms)
<a href="#">XCAMLITE4K8MPA</a>	Sony IMX334(C) 1/1.8"(7.68x4.32)	2.0x2.0	505mv with 1/30s 0.1mv with 1/30s	30@3840*2160(HDMI) 20@3840*2160(USB)	1x1	0.04~1000
<a href="#">XCAMLITE4K8MPB</a>	Sony IMX485(C) 1/1.2"(11.14x6.26)	2.9x2.9	2188mv with 1/30s 0.39mv with 1/30s	30@3840*2160(HDMI) 20@3840*2160(USB)	1x1	0.04~1000



Figure 2-1 Available Ports on the Back Panel of the Camera Body

## The XCAMLITE4K Series Camera Help Manual

Interface	Function Description
USB Mouse	Connect USB mouse for easy operation with embedded XCamView software;
USB Video	Connect PC or other host device to realize video image transmission;
HDMI	Comply with HDMI1.4 standard. 4K or 1080P format video output for standard monitor;
DC12V	Power adapter connection (12V/1A);
SD	Comply with SDIO3.0 standard and SD card could be inserted for video and images storage;
LED	LED status indicator;
ON/OFF	Power switch;
Video Output Interface	Function Description
HDMI Interface	Comply with HDMI1.4 standard; 30fps@4K or 30fps@1080P;
USB Video Interface	Connecting USB port of PC for video transfer ; MJPEG format video;
Other Function	Function Description
Video Saving	Video format: 8M(3840*2160) H264/H265 encoded MP4 file; Video saving frame rate: 30fps;
Image Capture	8M (3840*2160) JPEG/TIFF image in SD card ;
Measurement Saving	Measurement information saved in layer mode with image content; Measurement information is saved together with image content in burn in mode.
ISP Function	Exposure(Automatic / Manual Exposure) / Gain, White Balance(Manual / Automatic / ROI Mode), Sharpening, 3D Denoise, Saturation Adjustment, Contrast Adjustment, Brightness Adjustment, Gamma Adjustment, Color to Gray, 50HZ/60HZ Anti-flicker Function
Image Operations	Zoom In/Zoom Out, Mirror/Flip, Freeze, Cross Line, Overlay, Embedded Files Browser, Video Playback, Measurement Function
Embedded RTC(Optional)	To support accurate time on board
Restore Factory Settings	Restore camera parameters to its factory status
Multiple Language Support	English / Simplified Chinese / Traditional Chinese / Korean / Thai / French / German / Japanese / Italian / Russian
Software Environment under USB Video Output	
White Balance	Auto White Balance
Color Technique	Ultra-Fine™ Color Engine
Capture/Control SDK	Windows/Linux/macOS/Android Multiple Platform SDK(Native C/C++, C#/VB.NET, Python, Java, DirectShow, Twain, etc)
Recording System	Still Picture or Movie
Operating System	Microsoft® Windows® XP / Vista / 7 / 8 / 8.1 /10 /11(32 & 64 bit) OSx(Mac OS X) Linux
PC Requirements	CPU: Equal to Intel Core2 2.8GHz or Higher
	Memory: 4GB or More
	Display:19" or Larger
	CD-ROM
Operating Environment	
Operating Temperature (in Centidegree)	-10°~ 50°
Storage Temperature (in Centidegree)	-20°~ 60°
Operating Humidity	30~80%RH
Storage Humidity	10~60%RH

### 3 Dimension of XCAMLITE4K Series



图 3-1 Dimension of XCAMLITE4K Series

### 4 XCAMLITE4K Series Camera Packing Information(TBD)



Figure 4-1 The XCAMLITE4K Series Camera Packing Information

Standard Packing List		
A	Gift box : L:25.5cm W:17.0cm H:9.0cm (1pcs,1.48kg/ box)	
B	XCAMLITE4K series camera	
C	Power Adapter: Input: AC 100~240V 50Hz/60Hz, Output: DC 12V 1A American standard: Model: GS12U12-P11 12W/12V/1A; UL/CUL/BSMI/CB/FCC European standard: Model: GS12E12-P11 12W/12V/1A; TUV(GS)/CB/CE/ROHS EMI Standard: EN55022, EN61204-3, EN61000-3-2, -3, FCC Part 152 class B, BSMI CNS14338 EMS Standard: EN61000-4-2, 3, 4, 5, 6, 8, 11, EN61204-3, Class A Light Industry Standard	
D	USB Mouse	
E	HDMI cable	
F	USB2.0 A male to A male gold-plated connectors cable /2.0m	
G	CD (Driver & utilities software, Ø12cm)	
Optional Accessory		
H	SD card(16G or above; Speed: class 10)	
I	Adjustable lens adapter	C-Mount to Dia.23.2mm eyepiece tube (Please choose 1 of them for your microscope) 108001/AMA037 108002/AMA050 108003/AMA075
J	Fixed lens adapter	C-Mount to Dia.23.2mm eyepiece tube (Please choose 1 of them for your microscope) 108005/FMA037 108006/FMA050 108007/FMA075
<b>Note:</b> For I and J optional items, please specify your camera type(C-mount, microscope camera or telescope camera), ToupTek engineer will help you to determine the right microscope or telescope camera adapter for your application;		

<b>K</b>	108015(Dia.23.2mm to 30.0mm Ring)/Adapter rings for 30mm eyepiece tube	
<b>L</b>	108016(Dia.23.2mm to 30.5mm Ring)/ Adapter rings for 30.5mm eyepiece tube	
<b>M</b>	Calibration kit	106011/TS-M1(X=0.01mm/100Div.); 106012/TS-M2(X,Y=0.01mm/100Div.); 106013/TS-M7(X=0.01mm/100Div., 0.10mm/100Div.)

## 5 Software and App

The software or the APP can be downloaded from the following link:

Windows: <http://www.touptek.com/download/showdownload.php?lang=en&id=33>

Linux & macOS: <http://www.touptek.com/download/showdownload.php?lang=en&id=28>

## 6 XCAMLITE4K Series Camera Application Configurations

You can use the XCAMLITE4K series camera in two different ways. Each application requires different hardware environment.

### 6.1 Camera Working Standalone with Built-in XCamView Software



Figure 6-1 XCAMLITE4K Series Camera with the HDMI Displayer

For this application, apart from the microscope, the user only needs an [XCAMLITE4K](#) series camera, an HDMI display, an HDMI cable, an SD card, a USB mouse and a power adapter that come with the camera. The steps to start the camera are listed as below:

- Connect the camera to a HDMI display using the HDMI cable;
- Insert the supplied USB mouse to the camera's USB port;
- Insert the supplied SD card into the [XCAMLITE4K](#) series camera SD card slot;
- Connect power adapter to the camera the and switch it on.;
- Switch on the display and view the video in the [XCamView](#) software. Move the mouse to the left, top or bottom of the [XCamView](#) UI, different control panel or UI will pop up and users could operate with the mouse at ease.

### 6.2 Connecting camera to the PC with USB port

For Windows user (Windows XP (32bit), Windows 7/8/10 (32/64 bit) ), please use [ToupView](#).

For macOS and Linux user (macOS 10.10 or above or Linux distributions with kernel 2.6.27 or higher), please use [ToupLite](#).

The steps to start the camera are listed below:

- Install the [ToupView/ToupLite](#) on your PC;
- Connect power adapter to the camera the and switch it on. After starting the camera, plug one end of the USB cable into the USB 2.0 port of the [XCAMLITE4K](#) series camera, and plug the other end into the USB port of the PC;
- Open [ToupView/ToupLite](#) software. The [XCAMLITE4K](#) series camera will be recognized automatically by software. In [ToupView/ToupLite](#) software, select the corresponding [XCAMLITE4K](#) series camera by clicking the camera name in the camera list.

#### Note:

When the USB cable and the mouse are plugged into the camera at the same time, the USB cable is preferred and the mouse is not available; when the USB cable is unplugged, the mouse can be used normally.

## 7 Brief Introduction of XCAMLITE4K Series Camera's UI and Its Functions

### 7.1 XCamView UI



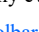





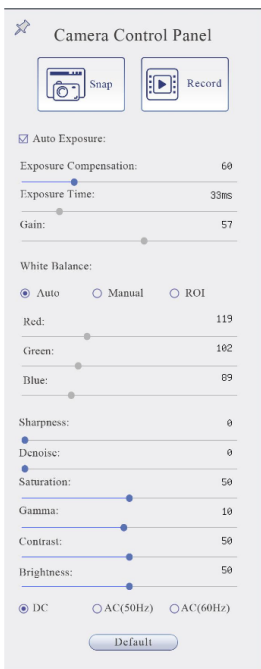
Figure 7-1 The XCAMLITE4K Series Camera Control GUI

Notes	
1	To show the <a href="#">Camera Control Panel</a> , move your mouse to the left of the video window. See <a href="#">Sec.7.2</a> for details.
2	Move the mouse cursor to the top of the video window, a <a href="#">Measurement Toolbar</a> will pop up for calibration and measurement operations. When user left-clicks the <a href="#">Float/Fixed</a> button  on the <a href="#">Measurement Toolbar</a> , the <a href="#">Measurement Toolbar</a> will be fixed. In this case the <a href="#">Camera Control Panel</a> will not pop up automatically even if users move mouse cursor to left side of the video window. Only when user left-clicks the  button on the <a href="#">Measurement Toolbar</a> to exit from measuring operations will they be able to do other operations on the <a href="#">Camera Control Panel</a> , or the <a href="#">Synthesis Camera Control Toolbar</a> . During the measuring operations, when a specific measuring object is selected, an <a href="#">Object Location &amp; Attributes Control Bar</a>  will appear for changing location and properties of the selected object. See <a href="#">Sec.7.4</a> for details.
3	When users move mouse cursor to the bottom of the video window, the <a href="#">Synthesis Camera Control Toolbar</a> will pop up automatically.  . See <a href="#">Sec.7.4</a> for details.



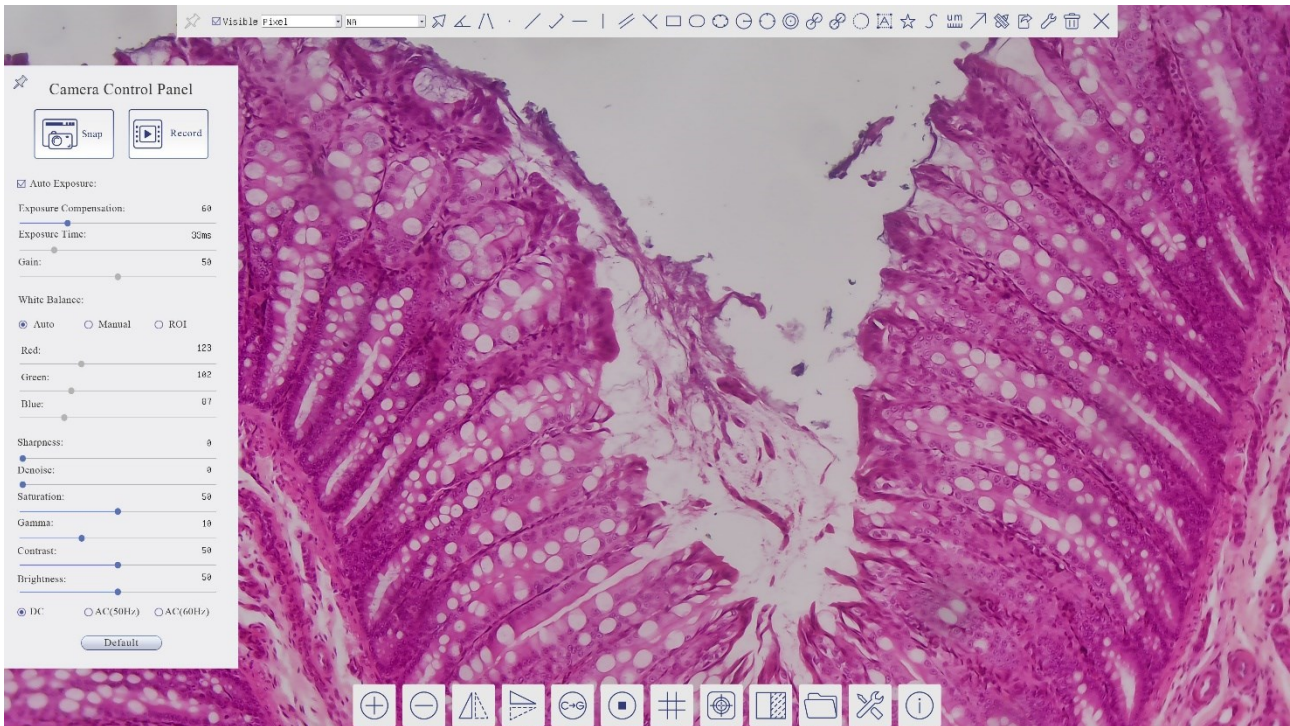




### 7.2 The Camera Control Panel on the Left Side of the Video Window

The **Camera Control Panel** controls the camera to achieve the best video or image quality according to the specific applications; It will pop up automatically when the mouse cursor is moved to the left side of the video window (in measurement status, the **Camera Control Panel** will not pop up. The **Camera Control Panel** will only pop up when the measurement operations are finished or terminated while user's cursor on the left edge of the video window). Left-clicking  button to achieve **Display/Auto Hide** switch of the **Camera Control Panel**.













Camera Control Panel	Function	Function Description
	<b>Snap</b>	<b>Capture</b> image and save it to the SD card
	<b>Record</b>	<b>Record</b> video and save it to the SD card
	<b>Auto Exposure</b>	When <b>Auto Exposure</b> is checked, the system will automatically adjust exposure time and gain according to the value of exposure compensation
	<b>Exposure Compensation</b>	Available when <b>Auto Exposure</b> is checked. Slide to left or right to adjust <b>Exposure Compensation</b> according to the current video brightness to achieve proper brightness value
	<b>Exposure Time</b>	Available when <b>Auto Exposure</b> is not checked. Slide to left or right to reduce or increase <b>Exposure Time</b> , adjusting brightness of the video
	<b>Gain</b>	Adjust <b>Gain</b> to reduce or increase brightness of video. The Noise will be reduced or increased accordingly
	<b>Auto White Balance</b>	<b>White Balance</b> adjustment according to the video continuously
	<b>Manual White Balance</b>	Adjust the <b>Red</b> or <b>Blue</b> slide bar to set the video <b>White Balance</b> .
	<b>ROI White Balance</b>	<b>White Balance</b> could be adjusted when the <b>ROI</b> region is changed according to content inside the <b>ROI</b> region.
	<b>Red</b>	Slide to left or right to decrease or increase the proportion of <b>Red</b> item in RGB on video
	<b>Green</b>	Slide to left or right to decrease or increase the proportion of <b>Green</b> item in RGB on video
	<b>Blue</b>	Slide to left or right to decrease or increase the proportion of <b>Blue</b> item in RGB on the video
	<b>Sharpness</b>	Adjust <b>Sharpness</b> level of the video
	<b>Denoise</b>	Slide left or right to <b>Denoise</b> the video
	<b>Saturation</b>	Adjust <b>Saturation</b> level of the video
	<b>Gamma</b>	Adjust <b>Gamma</b> level of the video. Slide to the right side to increase <b>Gamma</b> and to the left to decrease <b>Gamma</b> .
	<b>Contrast</b>	Adjust <b>Contrast</b> level of the video. Slide to the right side to increase <b>Contrast</b> and to the left to decrease <b>Contrast</b> .
	<b>DC</b>	For <b>DC</b> illumination, there will be no fluctuation in light source so no need for compensating light flickering
	<b>AC(50HZ)</b>	Check <b>AC(50HZ)</b> to eliminate flickering caused by 50Hz light source
	<b>AC(60HZ)</b>	Check <b>AC(60HZ)</b> to eliminate flickering caused by 60Hz light source
<b>Default</b>	Restore all the settings in the <b>Camera Control Panel</b> to default values	

### 7.3 The Measurement Toolbar on Top of the Video Window



The **Measurement Toolbar** will pop up when moving mouse cursor to any place near the upper edge of the video window. Here is the introduction of the various functions on the **Measurement Toolbar**:


Icon	Function
	<b>Float/ Fix</b> switch of the <b>Measurement Toolbar</b>
<input checked="" type="checkbox"/> Visible	Show / Hide <b>Measurement Objects</b>
Nanometer (nm)	Select the desired <b>Measurement Unit</b>
4X	Select <b>Magnification</b> for measurement after <b>Calibration</b>
	<b>Object Select</b>
	<b>Angle</b>
	<b>4 Points Angle</b>
	<b>Point</b>

## The XCAMLITE4K Series Camera Help Manual

	Arbitrary Line
	3 Points Line
	Horizontal Line
	Vertical Line
	3 Points Vertical Line
	Parallel
	Rectangle
	Ellipse
	5 Points Ellipse
	Circle
	3 Points Circle
	Annulus
	Two Circles and its Center Distance
	3 Points Two Circles and its Center Distance
	Arc
	Text
	Polygon
	Curve
	Scale Bar
	Arrow
	Execute <a href="#">Calibration</a> to determine the corresponding relation between magnification and resolution, which will establish the corresponding relationship between <a href="#">Measurement Unit</a> and the sensor pixel size. <a href="#">Calibration</a> needs to be done with the help of a micrometer. For detailed steps of carrying out <a href="#">Calibration</a> , please refer to <a href="#">ToupView</a> help manual.
	Export the Measurement information to CSV file(*.csv)
	Measurement Setup
	Delete all the measurement objects from
	Exit from <a href="#">Measurement</a> mode
	When the measurement ends, left-click on a single measuring object and the <a href="#">Object Location &amp; Properties Control Bar</a> will show up. User could move the object by dragging the object with the mouse. But more accurate movement could be done with the control bar. The icons on the control bar mean <a href="#">Move Left</a> , <a href="#">Move Right</a> , <a href="#">Move Up</a> , <a href="#">Move Down</a> , <a href="#">Color Adjustment</a> and <a href="#">Delete</a> .

### Note:

1) When user left-clicks [Display/Hide](#) button  on the [Measurement Toolbar](#), the [Measurement Toolbar](#) will be fixed. In this case the [Camera Control Panel](#) will not pop up automatically even if moving the mouse cursor to the left edge of the video window. Only when user left-click the  button on the [Measurement Toolbar](#) to exit from the measurement mode will they be able to doing other operations with the [Camera Control Panel](#) or the [Synthesis Camera Control Toolbar](#).

2) When a specific [Measurement Object](#) is selected during the measurement operation, the [Object Location & Attributes Control Bar](#)  will appear for changing the object location and properties of the selected objects.

7.4 Icons and Functions of the Synthesis Camera Control Toolbar at the Bottom of the Video Window



Icon	Function	Icon	Function
	Zoom In the video Window		Zoom Out the video Window
	Horizontal Flip		Vertical Flip
	Color/Gray		Video Freeze
	Display Cross Line		Overlay
	Compare Image with the current video		Browse Images and Videos in the SD Card
	Settings		Check the Version of XCamView

The Setting function is relatively more complicated than the other functions. Here are more details about it:

7.4.1 Setting>Measurement

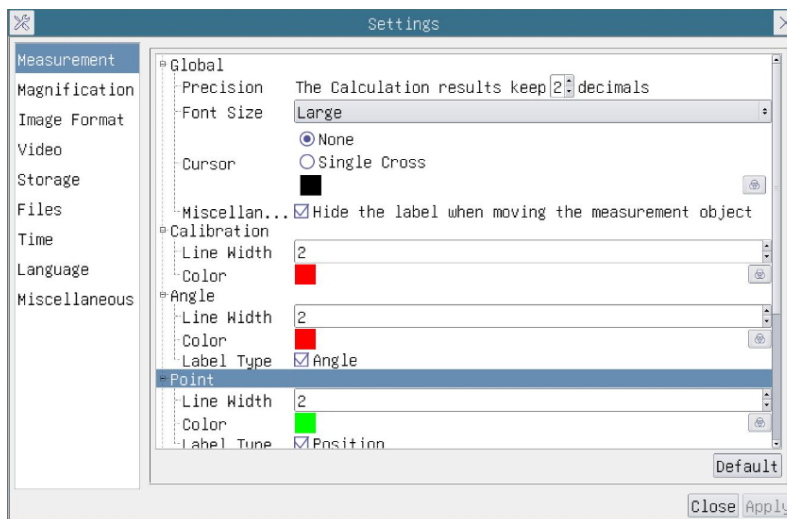


Figure 7-2 The Measurement Setup

Global	Precision	Used to set the number of digits after the decimal point of the measurement result
Calibration	Line Width	Used for defining width of the lines for Calibration;
	Color	Used for defining color of the lines for Calibration;
	EndPoint	Type: Used for defining shape of the Endpoint of lines for calibration: Null means no Endpoint, rectangle means rectangle type of Endpoint. It makes alignment more easily;
Point, Angle, Line, Horizontal Line, Vertical Line, Rectangle, Circle, Ellipse, Annulus, Two Circles, Polygon, Curve		
Left-click the  along with the Measurement command mentioned above will unfold the corresponding attribute settings to set the individual property of the Measurement Objects.		

7.4.2 Setting>Magnification

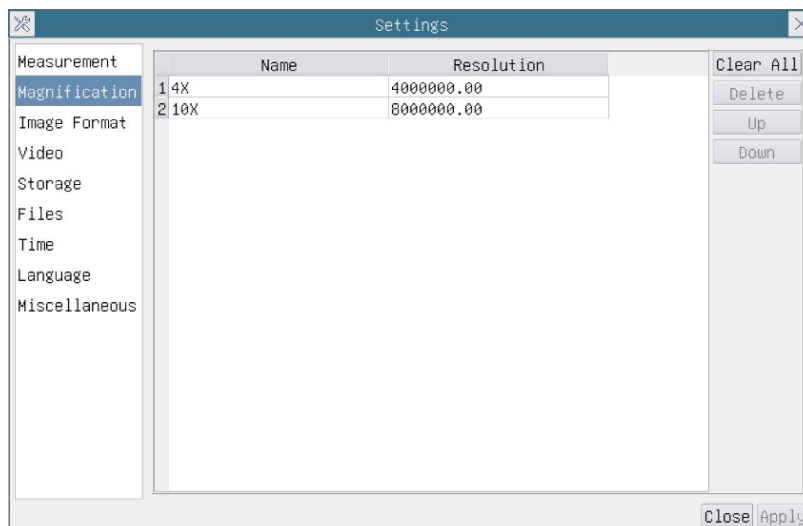


Figure 7-3 Comprehensive Magnification Calibration Settings Page

<b>Name</b>	The name of the <b>Magnification</b> , usually the <b>Magnification</b> of the objective of the microscope is used as the <b>Magnification</b> name when calibration, such as <b>4X</b> , <b>10X</b> , <b>40X</b> , <b>100X</b> , etc. Besides, other user-defined information could be added into the <b>Magnification</b> name too, for example, microscope model, operator name, etc.
<b>Resolution</b>	Pixels per meter. Image device like microscopes have high resolution value;
<b>Clear All</b>	Click the <b>Clear All</b> button will clear the calibrated <b>Magnification</b> ;
<b>Delete</b>	Click <b>Delete</b> to delete the selected <b>Magnification</b> ;
<b>Up</b>	Click <b>Move Up</b> to move up the selected <b>Magnification</b> ;
<b>Down</b>	Click <b>Move Down</b> to move the selected <b>Magnification</b> down;

7.4.3 Settings>Image Format

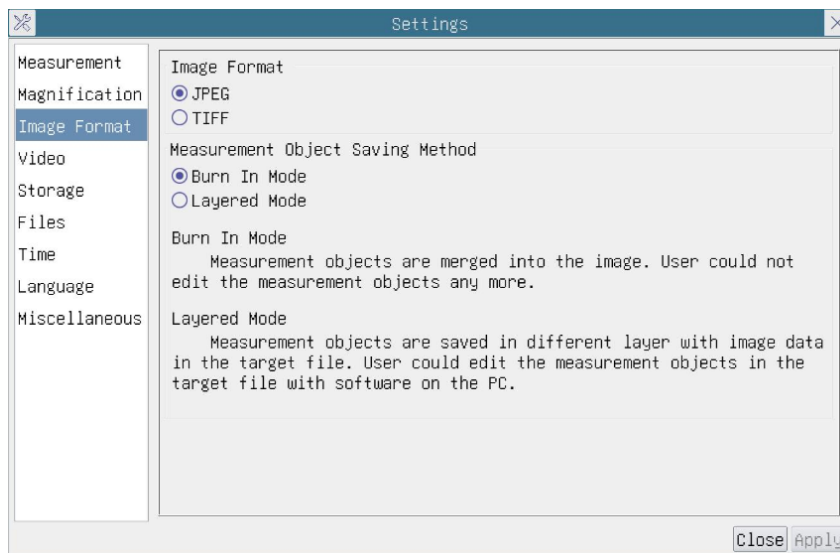


Figure 7-4 Comprehensive Image Format Settings Page

<b>Image Format</b>	<b>JPEG:</b> The extension of JPEG file can get very high compression rate and display very rich and vivid images by removing redundant images and color data. In other words, it can get better image quality with the least disk space. If <b>Measurement Objects</b> are available, the <b>Measurement Objects</b> will be burned into the image and the <b>Measurement Objects</b> cannot be edited. <b>TIFF:</b> Tag Image File Format(TIFF) is a flexible bitmap format that is mainly used to store images including photos and artistic images.
<b>Measurement Object Saving Method</b>	<b>Burn in Mode:</b> The <b>Measurement Objects</b> are merged into the current image. User could not edit the <b>Measurement Objects</b> anymore. This mode is not reversible. <b>Layered Mode:</b> The <b>Measurement Objects</b> are saved in different layer with current image data in the target file. User could edit the <b>Measurement Objects</b> in the target file with some software on the PC. This mode is reversible.

7.4.4 Settings>Video

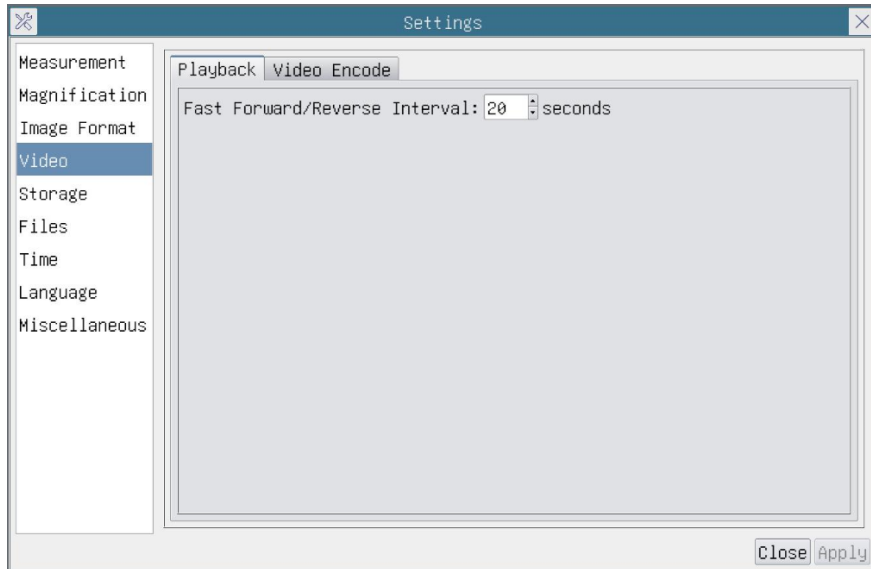


Figure 7-5 Comprehensive Setting of Video Settings Page-Playback

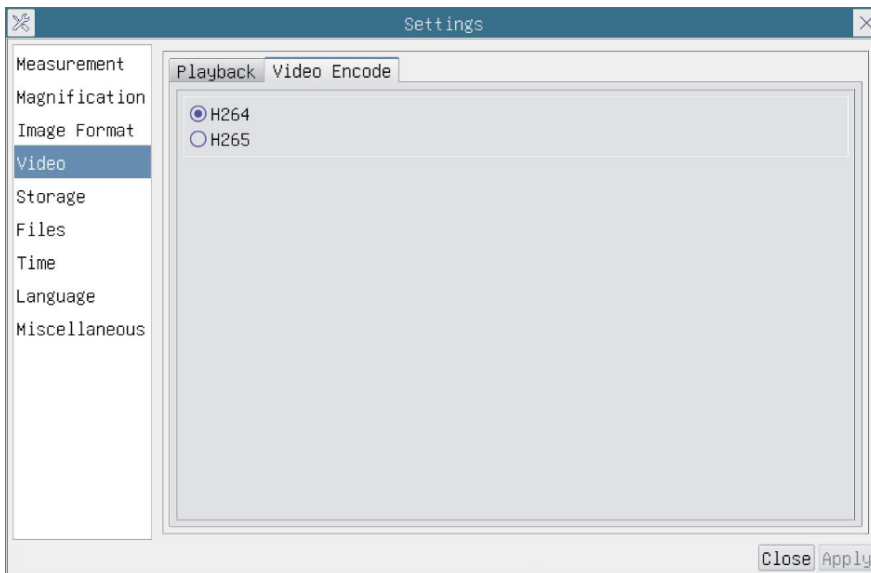


Figure 7-6 Comprehensive Setting of Video Settings Page-Video Encode

<a href="#">Fast Forward/Reverse Interval</a>	The time interval of the playback of video files.
<a href="#">Video Encode</a>	You can choose <a href="#">H264</a> or <a href="#">H265</a> encoding. <a href="#">H265</a> encoding can significantly reduce encoding bandwidth and save storage space under the same encoding quality;

7.4.5 Setting>Storage

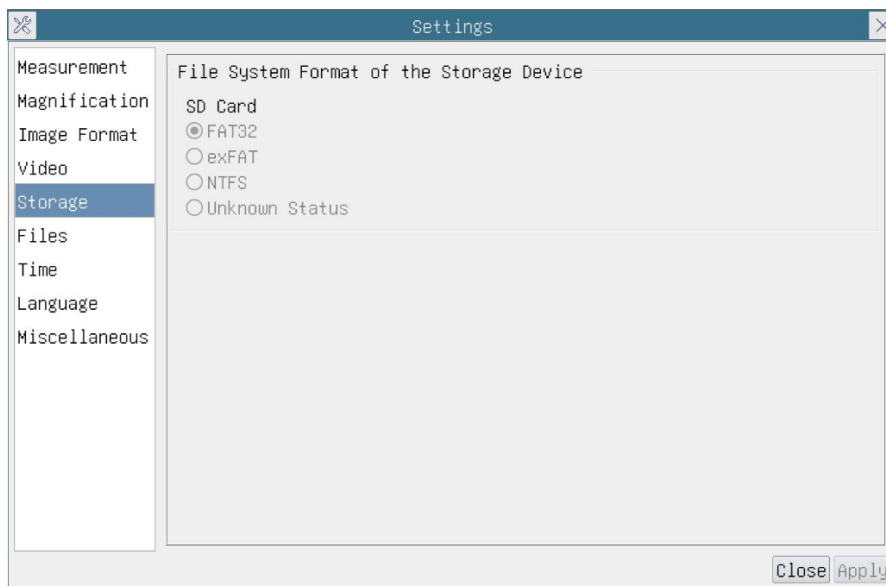


Figure 7-7 Comprehensive Setting of SD Card Setting Page

Storage Device	SD Card: SD Card is only supported as the storage device.
File System Format of the Storage Device	List the file system format of the current storage device FAT32: The file system of SD Card is FAT32. The maximum video file size of single file is 4G Bytes; exFAT: The file system of SD Card is exFAT. The maximum video file size of single file is 4G Bytes; NTFS: The file system of SD Card is NTFS. The maximum video file size of single file is 4G Bytes. Use PC to format the SD Card and switch between FAT32, exFat and NTFS. Unknown Status: SD Card not detected or the file system is not identified;

7.4.6 Setting>Files

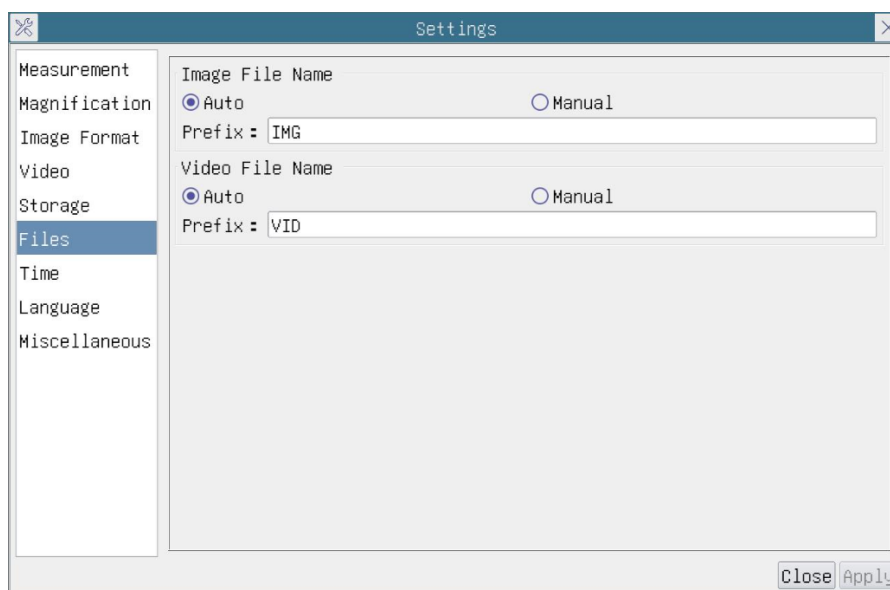


Figure 7-8 Comprehensive Setting of Files Settings Page

Image File Name	Auto: The image files will be saved automatically with the specified prefix. Manual: Users has to specify the file name before image saving.
Video File Name	Auto: The video file will be saved automatically with the specified prefix. Manual: Users has to specify the Video File Name before video recording.

### 7.4.7 Setting>Time

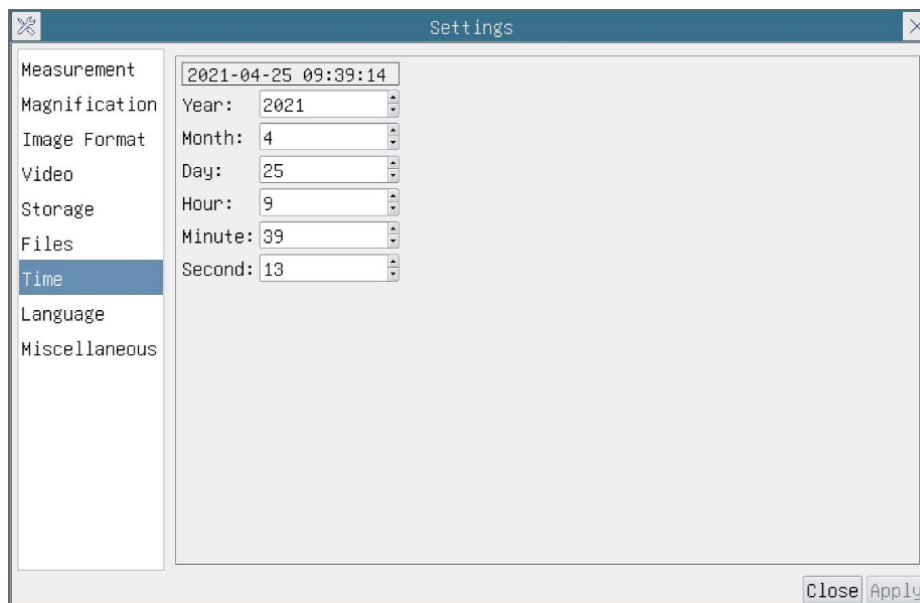


Figure 7-9 Time Setting

Time	User can set Year, Month, Day, Hour, Minute and Second in this page.
------	--

### 7.4.8 Setting>Language

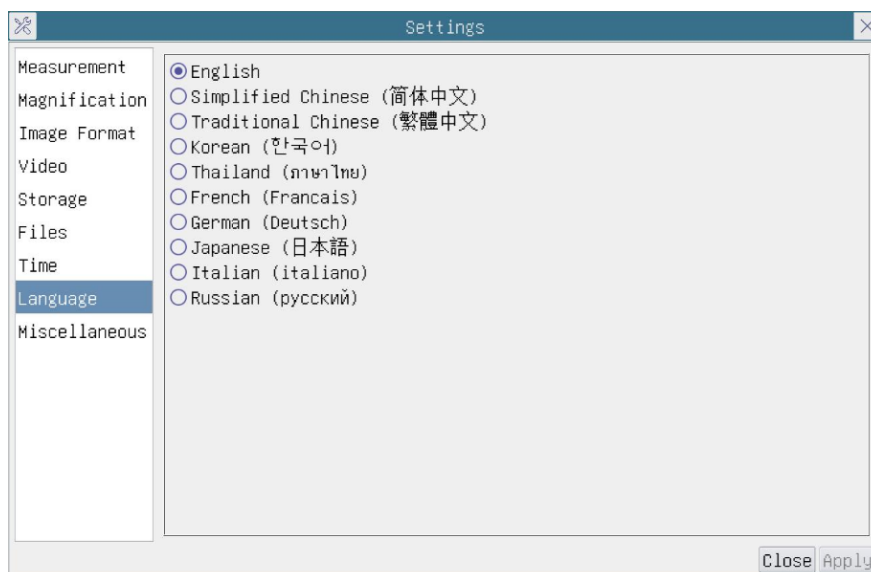


Figure 7-10 Comprehensive Setting of Language Selection Setting Page

English	Set language of the whole software into English;
Simplified Chinese	Set language of the whole software into Simplified Chinese;
Traditional Chinese	Set language of the whole software into Traditional Chinese;
Korean	Set language of the whole software into Korean;
Thailand	Set language of the whole software into Thailand;
French	Set language of the whole software into French;
German	Set language of the whole software into German;
Japanese	Set language of the whole software into Japanese;
Italian	Set language of the whole software into Italian;
Russian	Set language of the whole software into Russian;

7.4.9 Comprehensive Miscellaneous Settings Page

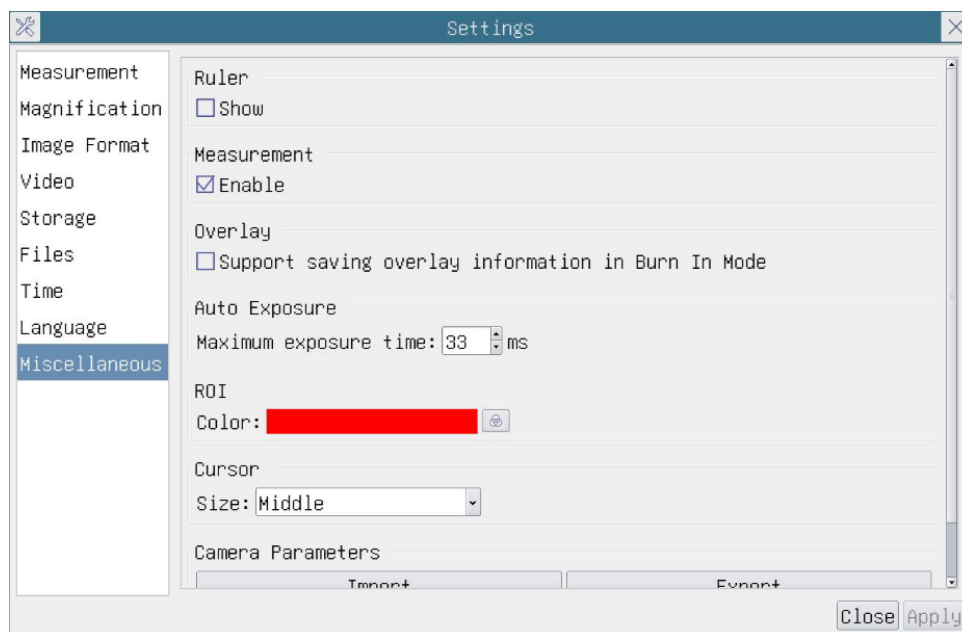


Figure 7-11 Comprehensive Miscellaneous Settings Page

<b>Ruler</b>	Select to display the <b>Ruler</b> in the video window, or not to display the <b>Ruler</b> ;
<b>Measurement</b>	Select to display the <b>Measurement</b> toolbar in the video window, otherwise, the <b>Measurement</b> toolbar will not be displayed;
<b>Overlay</b>	Select to support saving graphics <b>Overlay</b> information in fusion mode, and not to save graphics <b>Overlay</b> information in fusion mode;
<b>Auto Exposure</b>	The maximum exposure time during auto exposure process could be specified. Setting this item to a lower value could guarantee a faster frame rate during <b>Auto Exposure</b> ;
<b>ROI Color</b>	Choosing the <b>ROI</b> rectangle line color ;
<b>Camera Parameters Import</b>	Import the <b>Camera Parameters</b> from the SD card to use the previously exported <b>Camera Parameters</b> ;
<b>Camera Parameters Export</b>	Export the <b>Camera Parameters</b> to the SD card to use the previously exported <b>Camera Parameters</b> ;
<b>Reset to factory defaults</b>	Restore camera parameters to its factory status;



## 8 Sample Photos Captured with XCAMLITE4K Series Camera

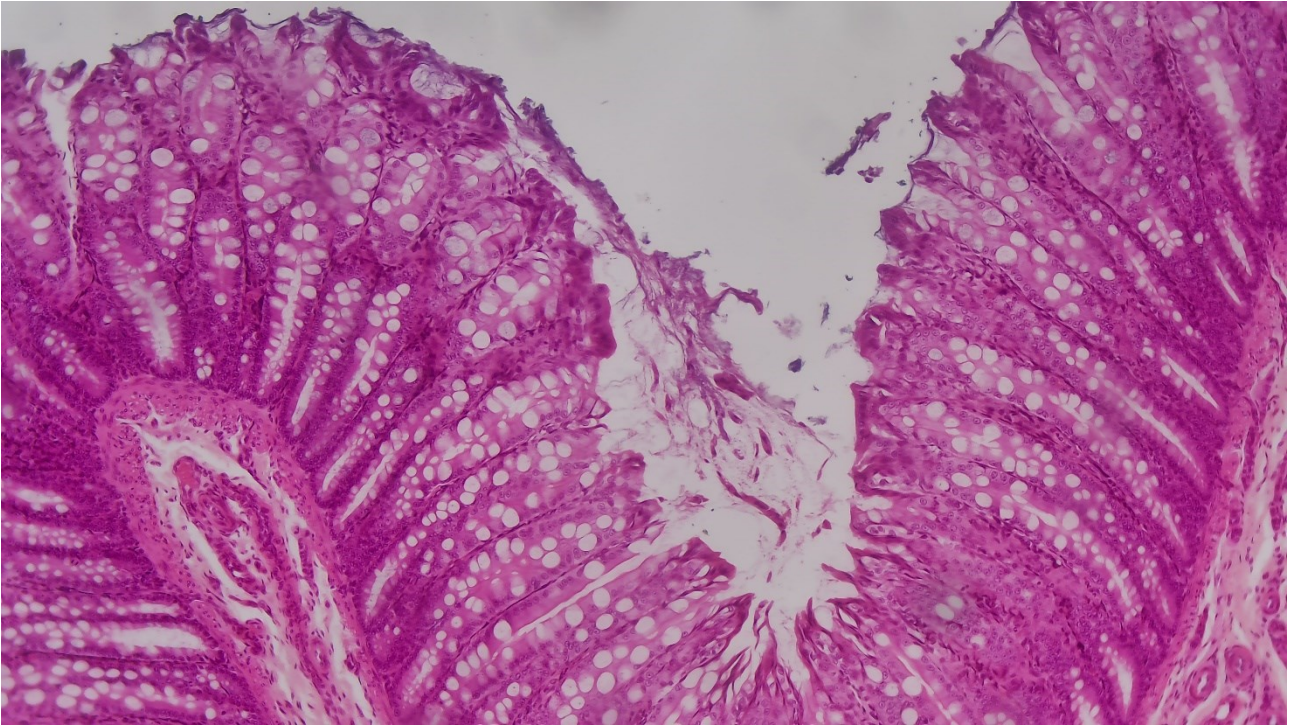


Figure 8-1 Large Intestine Captured with XCAMLITE4K8MPA

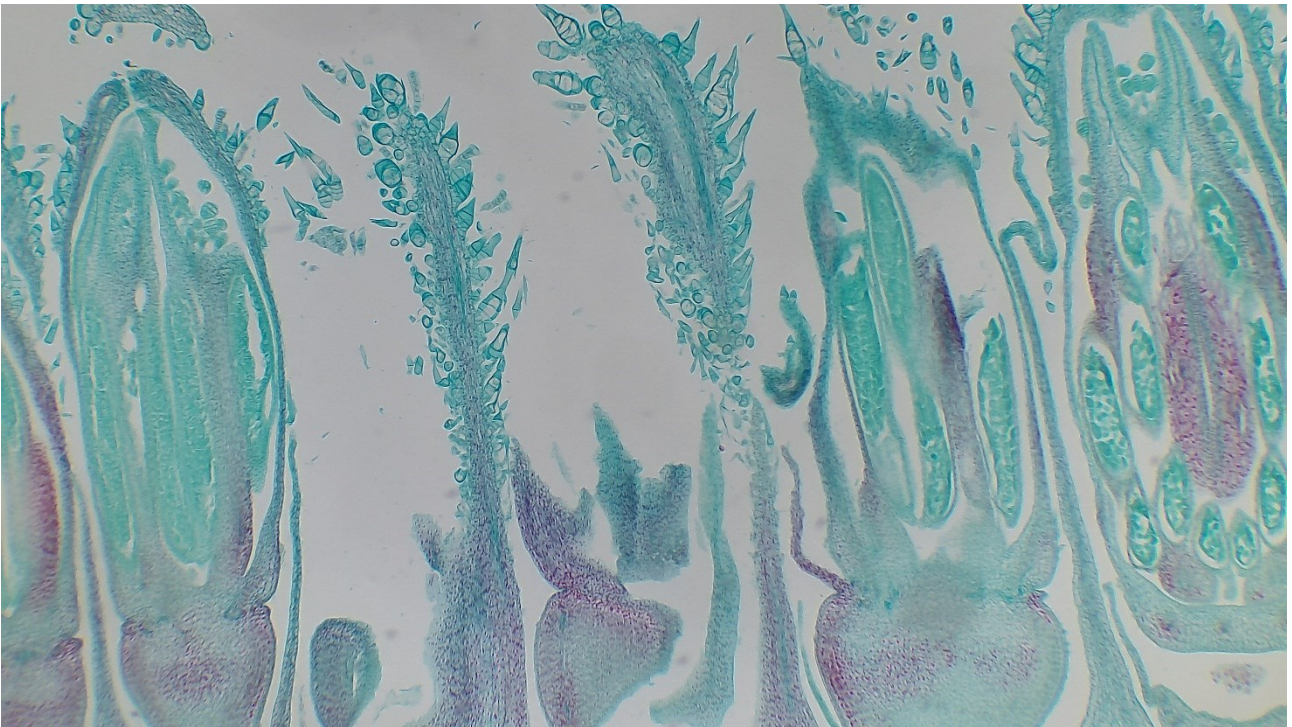


Figure 8-2 Sunflower Flower Captured with XCAMLITE4K8MPA

## 9 Contacting Customer Service

Please contact your local distributor if you have any questions about the product.